



Northern Lake Service Inc.

ENVIRONMENTAL ANALYTICAL LABORATORY

400 N Lake Ave – Crandon, WI 54520 | 2420 N Grandview Blvd – Waukesha, WI 53188

P: 800-278-1254 F: 715-478-3060 E: clientservices@nlslab.com

VOC, PVOC & GRO Solids/Soils Sample Collection

Contents of Sampling Kit:

- 2 x 2 oz (60mL) tared glass jars
- 2 x 40mL vials of methanol
- Methanol Trip Blank – DO NOT OPEN

Sample Collection:

1. Per method, weigh out approximately 25 grams of soil and deposit into a tared glass jar.
2. Pour one vial of methanol over the soil inside the tared glass jar.
3. Repeat with second set.
4. Upon successfully depositing the soil sample into the tared glass jar and adding the methanol, brush away any soil particles from the threads of the glass jar and the jar lid.