



Northern Lake Service Inc.

ENVIRONMENTAL ANALYTICAL LABORATORY

400 N Lake Ave – Crandon, WI 54520 | 2420 N Grandview Blvd – Waukesha, WI 53188

P: 800-278-1254 F: 715-478-3060 E: clientservices@nlslab.com

VOC, PVOC & GRO Aqueous Sample Collection

Contents of Sampling Kit:

- 2 x 40mL clear glass vials preserved with dilute HCL
- Trip Blank – DO NOT OPEN

Sample Collection:

1. Slowly fill each vial so that the water domes up just before it overflows. Avoid turbulence.
2. Replace the cap gently. When vials are inverted, there should be no visible air pockets or bubbles in the vials. (This condition is known as zero headspace)
3. If bubble(s) are present, uncap the vial and add a little more sample.