



# Northern Lake Service Inc.

## ENVIRONMENTAL ANALYTICAL LABORATORY

### **Fluoride Test Kit - \$31.50**

#### **Kit includes:**

- 1 x 125 mL Plastic Bottle

#### **General Sampling Instructions:**

**READ ALL INSTRUCTIONS BEFORE PROCEEDING WITH SAMPLE COLLECTION**

**\*Sampling video available on our website at [www.nlslab.com](http://www.nlslab.com) under support material**

Collect water that is representative of the way you actually consume your water.

1. Collect water from selected location, avoiding new plumbing if possible.
2. Remove aeration screen from faucet if one is present.
3. Run water at maximum flow for at least 2 minutes.
4. Fill the container.

#### **Follow the return shipping instructions to ensure kit is shipped back to the lab properly:**

- Fill out the required information below, and label your sample bottles with date/time of collection.
- Samples are accepted 7:00 am to 4:00 pm Monday – Friday at the Crandon lab and 8:00 am to 4:00 pm Monday – Friday at the Waukesha lab. If you are shipping your kit to NLS, samples **MUST** arrive at the laboratory no longer than 28 days from the time of collection.
- NLS receives deliveries from UPS, Fed Ex, SpeeDee Delivery, USPS, and other private courier services. Use any of these courier services to return sample to NLS.
- If you have any questions regarding this sampling kit, please call our Client Service Department at 800-278-1254

# Fluoride Test Kit - \$31.50

## Return your sample as soon as possible to one of our two locations:

Northern Lake Service  
400 N Lake Ave  
Crandon, WI 54520

**OR**

Northern Lake Service  
2420 N Grandview Blvd  
Waukesha, WI 53188

## Please provide the following information:

Name: \_\_\_\_\_

Address: \_\_\_\_\_

City/State/Zip: \_\_\_\_\_

Phone: \_\_\_\_\_ Email: \_\_\_\_\_

Sample Collection Date: \_\_\_\_\_ Sample Collection Time: \_\_\_\_\_ AM/PM

Sample Collection Location (ex. Pressure Tank) \_\_\_\_\_

Sample Collected By: \_\_\_\_\_

Report by email only  Report by mail only  Report by email and mail